

Exporting Images

In OMERO you are able to export any image in OME-tiff, save it as a JPEG, or if it was archived at import, download it in the original image format.

1 Click on the **Saving Options** icon in the toolbar.



2 If the image was archived in its original file format select **Download** from the dropdown menu.

			Download Image: Comparison of Co
	Download	Download. Eports sOMC-TFF Image 1D: 73 F-TEE_19_B3D_D3D.dv collected with critical illumination	Save in: Dourhoods
		Download	Network Files of type: Al Files

3 Select **Export as OME-TIFF** from the dropdown menu to export in that format.







4 Select **Save as JPEG** to export as a JPEG, PNG or TIFF.

.	Save as JPEG	Collected with critical illumination	→	Save As Set of where to save locally the mages as 3PEG. Save As Save As Save As Prite Prit
		Save		Corputer Computer Reder name: *Documents[1 CMET/restrupTrest Data 2012/Data from Jab Network Ples of type: PEG (*Jong, *Jog, *Jog) ftum#able Save

5 Scripts can be used to output in a number of formats.



